



**Newport.**

**ORIEL PRODUCT LINE**

150 Long Beach Boulevard  
Stratford, CT 06615

Phone: (203) 377-8282  
(800) 714-5393

Fax (203) 378-2457

E-MAIL: [oriel.sales@newport.com](mailto:oriel.sales@newport.com)

**MS260i™**

**IMAGING 1/4 M SPECTROGRAPH**

**MODELS 74050, 74055, 74085, 74086**

**USER MANUAL**

Please read these instructions completely before operating this equipment. The specification and operating instructions apply only to the model(s) covered by this manual. If there are any questions or problems regarding the use of this equipment, please contact Newport or the representative from whom this equipment was purchased.

**Rev: 04-24-09**

## TABLE OF CONTENTS

I. INTRODUCTION.....	3
I.1 OPTICAL CONFIGURATION.....	3
I.2 DESIGNED FOR MULTI-TRACK SPECTROSCOPY .....	3
I.3 NEGLIGIBLE STRAY LIGHT .....	3
I.4 EXCELLENT SPECTRAL RESOLUTION .....	3
I.5 WAVELENGTH DRIVE .....	3
II. UNPACKING AND SETTING UP .....	4
II.1 SOFTWARE INSTALLATION .....	5
II.2 USB INSTALLATION FOR THE 74125.....	<b>ERROR! BOOKMARK NOT DEFINED.</b>
II.3 UTILITY PROGRAM INSTALLATION.....	6
III. SAFETY .....	8
IV. 15 MINUTE BASICS.....	9
IV.1 GETTING STARTED WITH THE HAND CONTROLLER .....	9
IV.2 GETTING STARTED WITH IEEE-488.....	12
IV.3 GETTING STARTED WITH RS-232 COMMUNICATIONS.....	13
V. FUNCTIONS AND CONTROLS .....	15
V.1 COMMAND AND QUERY SYNTAX .....	15
V.2 HANDSHAKING MODE.....	16
V.3 BRIEF COMMAND REFERENCE ORGANIZED ALPHABETICALLY .....	17
V.4 DETAILED COMMAND REFERENCE ORGANIZED BY FUNCTION.....	18
V.5 ERROR CODES WITH DESCRIPTIONS .....	23
VI. MS260i™ UTILITY PROGRAM .....	24
VI.1 SOURCE CODE .....	25
VI.2 IEEE-488 VERSION .....	25
VI.3 RS-232 VERSION .....	26
VI.4 USB VERSION 74085/74086 .....	26
VII. OPTIONAL ACCESSORIES .....	27
VII.1 MICROMETER DRIVEN SLIT .....	27
VII.2 FILTER WHEEL.....	27
VII.3 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION.....	27
VII.4 MEMORY LOSS (LARGE WAVELENGTH CALIBRATION ERRORS) .....	31
VIII. DECLARATION OF CONFORMITY .....	32
IX. WARRANTY AND RETURNS .....	33

## I. INTRODUCTION

---

This new spectrograph is in answer to your requests for an economical, fully automated, multi-grating, ¼ m imaging spectrograph. For those familiar with our popular MS257™ ¼ m Spectrograph, the MS260i™ meets many of the performance standards in a simplified design, and is more economical.

The MS260i™ offers the best performance to price ratio of any spectrograph on the market.

The MS260i™ is an F/3.9 instrument with a single flat, 28 x 10 mm output port for multichannel detectors. Stray light is negligible, throughput is high and spatial and spectral resolution are impressive. Three communication interfaces are available via two versions of the board; one with RS232 and IEEE488 and one with USB only. Control via optional Hand Controller is available.

This manual applies to the:

- 74050 MS260** MS260 (single output port) which has **RS232 and GPIB** communications interface
- 74085 MS260** MS260 (single output port) which has a **USB** communications interface
- 74055 MS260** MS260 (dual output port) which has **RS232 and GPIB** communications interface
- 74086 MS260** MS260 (dual output port) which has a **USB** communications interface

### I.1 OPTICAL CONFIGURATION

The MS260i™ features an asymmetrical in-plane Czerny-Turner optical configuration. The input F/# is F/3.9, and the input and exit focal lengths are unequal 220 and 257 mm. A set of computer designed toroidal mirrors produce accurate images of the input slit in the flat output plane.

### I.2 DESIGNED FOR MULTI-TRACK SPECTROSCOPY

Multi-track Spectroscopy is the simultaneous recording of multiple spectra. You can't find a better Multi-track Spectroscopy instrument combination than Oriel's Imaging CCD Detectors and MS260i™. One of the factors limiting the number of spectra, which can be recorded simultaneously, is the spectrograph's magnification. MS260i™ uses two toroidal mirrors to achieve optimum imaging. These mirrors have different horizontal and vertical radii. The result is a 1.1x horizontal and 1.6x vertical magnification.

### I.3 NEGLIGIBLE STRAY LIGHT

One of the design criteria for MS260i™ was to eliminate light scatter inside the instrument, one of the main contributors to stray light. It is important that diffracted light not be allowed to reflect from the mirrors, or from the face of a focal plane detector, back onto the grating. This light could then be re-diffracted and cause ghost images or stray light at the detector. The wide open optical layout and exaggerated tilt of the exit focal plane of the MS260i™ ensure that there is zero re-entrant spectra.

### I.4 EXCELLENT SPECTRAL RESOLUTION

Resolution of 0.25 nm, with a 0.25 µm slit and a 1200 l/mm grating, is possible with the MS260i™! This compares to 0.4 nm with a 25 µm slit for the MS127i™, 1/8 m Imaging Spectrograph.

### I.5 WAVELENGTH DRIVE

The MS260i™ uses a high speed stepping motor drive for rapid wavelength changes. This same drive is used to change the grating; a grating change takes only 10 seconds.

## II. UNPACKING AND SETTING UP

---

Carefully unpack the MS260i™ Spectrograph and its accessories. The packaging is molded to your instrument and therefore provides optimal protection. You should save the packaging in case you ever need to transport or ship this instrument.

Your MS260i™ comes completely assembled with up to three gratings installed. Your fixed slit housings, or micrometer driven slits, are mounted at our factory. The instrument is closed and the optical ports are covered to prevent contamination of the optics during shipment. Included with the MS260i™ you should also find:

1. this manual
2. a tabletop power supply (universal input)
3. a line cord for your country
4. any additional gratings you ordered
5. a CD-ROM containing your MS260i™ software

Since the MS260i™ is aligned and calibrated at our factory, there is no need to remove the top. You can get started immediately by connecting the 5 pin DIN connector of the power supply to the MS260i™ back panel, connect the line cord to both the power supply and the wall, and then turn on the MS260i™ power switch. You will hear the monochromator drive to its home position, then any accessories like Filter Wheel and/or Motorized Slits will home, finally the wavelength drive will position to the previous grating and wavelength position.

The external BNC on the chassis is to facilitate external TTL control of the Shutter. A TTL low (or short) will activate (Open) the Shutter. Control can be overridden by the shutter commands from a computer or the shutter key on the hand controller.

## II.1 SOFTWARE INSTALLATION

Newport provides an elementary software package with your Cornerstone™ MS260. To load and run the software you must have a PC compatible computer (with CDROM drive) running Microsoft Windows XP SP1 or above. For the 74085/74086 a USB2 port is required. **To successfully complete the install of the MS260 USB driver or the Utility software you must have Administrator privileges on the PC which you are performing the install.**

The CD contains the installation for the Utility program, LabView VIs and the USB drivers. The Utility Program provides easy access to almost all Cornerstone™ functions and is discussed in Section VI CORNERSTONE™ UTILITY PROGRAM. Additionally, you can use the LabView VIs when writing your own LabView software for Cornerstone™ (National Instruments LabView 8.5 required). The source code for both Utility program is included with the 74100 VIs to be used as an example.

There are three versions of the basic Cornerstone™ software. The first version is for IEEE-488 communications. This version requires that you have a National Instruments IEEE-488 controller card and the corresponding software drivers installed in your PC. The second version is for RS-232 communications. There is also a USB version for the 74125 version of the MS260, the USB driver must be installed for that version to function. In all cases, the software looks the same, and has the same functionality.

See the Table I below for where to find the Utility install program, the VIs and the USB drivers.

**Table I(Cornerstone Utility CD 74050/74055 74085/74086 relevant directories)**

	Directory on Utility CD	Sub-directory containing Install Program (Setup.exe) for the Utility application	Sub-directory containing LabView VIs
MS260 using RS232 Single output port	\740500RS232\	\74050RS232\Utility\Installer	\74050RS232\
MS260 using GPIB Single output port	\74050GPIB\	\74050GPIB\Utility\Installer	\74050GPIB\
74085 CS260 USB Single output port	\74085\	\74085\ Utility\Installer	\74085\
MS260 using RS232 Dual output port	\74055RS232\	\74055RS232\Utility\Installer	\74055RS232\
MS260 using GPIB Dual port output port	\74055GPIB\	\74055GPIB\ Utility\Installer	\74055GPIB\
74086 CS260 USB Dual output port	74086\	\74086\ Utility\Installer	\74086\
74085/74086 USB drivers and install file	\USB DRIVERS	--	--

## II.2 UTILITY PROGRAM INSTALLATION

**IMPORTANT:** Do not connect any USB Cornerstone™ to the computer unless the USB driver has been installed, or until prompted to do so.

**Note:** If the Utility software will not be installed with USB Model 74085/74086, please skip to §II.3 , p. 7.

1. Put the Cornerstone™ CD-ROM into your computer drive.
2. Select Run... from the Windows Start Menu. Then click on the Browse...button. Now use the controls in the Browse dialog box to select the Cornerstone™ CD, browse to the folder for your instrument as shown in Table I above. Inside the folder you will find "Setup.exe". Double-click on this program and then press the OK button in the Run dialog box.
3. The Installation program will appear, follow the prompts to install the Utility software on you PC. The default directories and Start Menu groups will work on most machines and should be used if possible. This simplifies the installation as you only need to click the "Next" button a few times.
4. Another install program will run and you should be able to use the default values again. Simply click "Finish" to complete the installation.
5. *If using USB Model 74085/74086, follow the connection procedure in §II.3 , p. 7 starting at Step #6.*
6. You are now ready to work with the Cornerstone™. Please review the Safety section in detail before proceeding to the 15 Minute Basics.

**<< IMPORTANT NOTE >>**

For the software to send proper numeric format to the instrument, Windows Regional Settings in Control Panel must be set to English (United States).

7. Put the Cornerstone™ CD-ROM into your computer drive.
8. Select Run... from the Windows Start Menu. Then click on the Browse...button. Now use the controls in the Browse dialog box to select the Cornerstone™ CD, browse to the folder for your instrument as shown in Table I above. Inside the folder you will find "Setup.exe". Double-click on this program and then press the OK button in the Run dialog box.
9. The Installation program will appear, follow the prompts to install the Utility software on you PC. The default directories and Start Menu groups will work on most machines and should be used if possible. This simplifies the installation, as you only need to click the "Next" button a few times.
10. Another install program will run and you should be able to use the default values again. Simply click "Finish" to complete the installation. You are now ready to work with the Cornerstone™. Please review the Safety section in detail before proceeding to the 15 Minute Basics.

### II.3 DRIVER INSTALLATION FOR USB MODEL 74085/74086

**IMPORTANT:** *Do not connect any USB Cornerstone™ to the computer unless the USB driver has been installed, or until prompted to do so.*

**Note:** If the monochromator(s) will be used with the optional TRACQ Basic™ software, follow its installation procedure now and disregard the remainder of this section.

Perform the following steps to install the USB driver for use with OEM or any other applications:

1. Log into a Windows user account *with Administrator privileges.*
2. Insert the Cornerstone™ CD-ROM into a CD or DVD drive and locate the USB drivers folder from Table I (§**Error! Reference source not found.**, p. **Error! Bookmark not defined.**).
3. Right-click on the file “oriel\_cs\_usb.inf” and select the *Install* option from the pop-up context menu.

**Note:** If filename extensions are not displayed, two files will appear as “oriel\_cs\_usb”. If the menu does not display the *Install* option for one file, use the other one.

4. If any warning dialogs appear stating that the driver is “not signed” or “has not passed Windows Logo testing”, click “Continue Anyway” to proceed.
5. Driver installation will complete silently if successful.

**Note:** If an “Installation failed” dialog appears, the logged-in Windows user account may not have Administrator privileges required for installation. *Please consult with your system manager and repeat from Step #1 before contacting Newport/Oriel Customer Service.*

6. Power-on one Cornerstone™ monochromator, then connect its USB cable to the computer. *Do not connect any additional devices yet.*
7. When the “Found New Hardware” wizard appears, select the option “Install the software automatically (Recommended)”, then click the “Next >” button. (It is not necessary to have the CD-ROM in the drive at this point.)
8. If any warning dialogs appear stating that the driver is “not signed” or “has not passed Windows Logo testing”, click “Continue Anyway” to proceed.
9. After the wizard completes, click the “Finish” button.
10. Repeat from Step #6 for any additional monochromators.

**Note:** If a Cornerstone™ is connected to a different USB port on the same computer at a later time, it may be necessary to repeat installation Steps #7 through #9.

### **III. SAFETY**

---

#### **Electrical**

The MS260i™ requires only DC voltages for operation. These voltages are provided by an external power supply. This universal input power supply accepts 110/220 VAC for operation and has no user serviceable parts. Do not attempt to open the external power supply. Do not attempt to work in the MS260i™'s electrical compartment without first disconnecting the power cord, since you may contact live voltage areas inside the compartment even with the power switch in the "off" position.

MS260i™ has an internal microprocessor and should be installed with appropriate surge/EMI/RFI protection on all power lines. Dedicated power lines or line isolation may be required for some extremely noisy sites. The circuits used in MS260i™ are extremely sensitive to static electricity and radiated electromagnetic fields. Operation of MS260i™ close to intense pulsed sources (lasers, xenon strobes or arc lamps, etc.) may compromise performance, if shielding is inadequate, and may cause permanent damage to the microprocessor. Note: This instrument conforms to European CE standards for both safety and EMC.

#### **Mechanical**

Avoid dropping, sudden shocks, or rough handling of the spectrograph since this may cause the system to lose its calibration and may destroy the high precision drive components or optics.

Do not use more than finger force in tightening down the grating mounts, since this may cause damage to the drive assembly.

#### **Optical**

Do not touch any optical surfaces since this is likely to cause irreparable damage. Do not attempt to clean any optical surface except by blowing off dust or lint particles with a stream of dry clean air or nitrogen.



## IV. 15 MINUTE BASICS

---

### IV.1 GETTING STARTED WITH THE HAND CONTROLLER

The 74009 Hand Controller is a fast and easy way to begin immediate communications with the MS260i™. Although the Hand Controller does not provide access to all the functions of this instrument, it does provide a simple interface to the most commonly used parameters and controls.

#### Physical Connections

Please note: None of the MS260i™ accessories are designed to be connected or disconnected with power on. Always turn off the instrument using the power switch on the back panel before you connect or remove the filter wheel or any communications cable (including the Hand Controller).

To get started with 74009 Hand Controller, simply connect the black communications cable by snapping the RJ-11 connector into the receptacle on top of the controller and snapping the other end into the matching connector on the back panel of your MS260i™.

#### Color Coding

The functional keys on the Hand Controller are divided into various color-coded groups as follows:

Red key at top left	“Abort” wavelength motion in progress, also clears pending commands.
Blue keys at the top	Control wavelength motion.
Red frame in row #2	Pertain to constant rate scanning.
Orange frame in row #2	Grating position and parameters.
Green frame on left side	Control accessories including filter, shutter, output port (when available).
Blue frame on left side	Toggle local/remote communications and sets operating units (shifted).

#### Activating the Hand Controller

With the Hand Controller connected, turn on power to the MS260i™. You will see a horizontal cursor in the upper left corner of the display. Press the LOCAL key to activate communications through the Hand Controller. Please note that if the Hand Controller is active (the display is lit and information fills the screen), you cannot command the MS260i™ via either RS-232, IEEE-488 or USB. Press the LOCAL key again to deactivate the Hand Controller, allowing PC communications.

#### Yellow “Shifted” Keys

Some keys are divided into top and bottom halves. The function for the yellow top half is called a “shifted” function. To activate these commands press the SHIFT key in the bottom left corner and then press the function key.

#### Entering Parameters- the Numeric Keypad

Many of the MS260i™ commands require a numeric parameter. First press the function key for your command. The bottom line of the LCD display will then show an abbreviated form of the command and the current parameter. There is no need to erase this number, simply typing the first key for your desired parameter will overwrite the current value. To execute the command press the ENTER key in the bottom right corner.

While entering parameters, if you mistype one digit press SHIFT and then DEL (above the decimal point in the numeric keypad); this will move backward one display position and delete that character.

---

## Key Reference

(From Top of the Controller to the Bottom, Left to Right)

- ABORT** If you accidentally press a function key or enter a bad parameter you may press the ABORT key to clear that command before it executes. The ABORT key has no effect on most functions once the command is executed (by pressing ENTER). However, wavelength motion will stop immediately upon ABORT and the display will update to show the wavelength where motion was halted.
- STEP DOWN** Press this key once to move the MS260i™ to shorter wavelengths by the smallest possible increment. Holding this key causes rapid single stepping toward lower wavelengths. Press SHIFT and then STEP DOWN, you can use the keypad to define a number of steps to move at once. Press ENTER to execute this larger move.
- STEP UP** Press this key once to move the MS260i™ to higher wavelengths by the smallest possible increment. Holding this key causes rapid single stepping toward higher wavelengths. Press SHIFT and then STEP UP, you can use the keypad to define a number of steps to move at once. Press ENTER to execute this larger move.
- GO WAVE** This key induces wavelength motion to a specified position. Press GO WAVE, then use the keypad to enter the desired wavelength, and finally press ENTER to execute the move.
- GO SCAN** Pertains to constant rate motion. This function is not available in the current version.
- SCAN  $\lambda$**  Pertains to constant rate motion. This function is not available in the current version.
- SHIFT SPEED** Pertains to constant rate motion. This function is not available in the current version.
- GRAT** Press this key to move to the other grating position. Press GRAT, then press either 1, 2 or 3 on the numeric keypad, and finally press ENTER to execute the change. The display will update to show the correct parameters for the new grating.
- SHIFT LABEL** This key (located directly above GRAT) provides access to the grating label. The label is an eight character alpha-numeric field that is used for information only. Newport sets the grating blaze wavelength as the label during calibration. However, you can use the keypad to type any number, then press ENTER to accept that new value. The Hand Controller display only shows the first four digits of the label, and you can only enter numbers.
- LINES** Press this key to change the lines/mm of the grating. This value is used for wavelength calculations and will affect your calibration. This parameter should be changed only when adding a new grating.
- SHIFT CALIB** This button is used to change the wavelength calibration of the current grating. Use the GO WAVE, STEP DOWN, and STEP UP keys to move to a known wavelength (often the intensity peak for a spectral calibration lamp). If the wavelength shown in the top row of the display does not match the known value, press SHIFT CALIB and use the numeric keypad to enter the known wavelength. Then press ENTER to change the calibration. The current grating is thus calibrated over its entire range.
- SHUTTER** Alternately open and close the shutter by pressing this button. The shutter position is shown in the display as "OPN" for open and "CLS" for closed.

- SHIFT PORT-** Use this key to alternate between the two output ports. This feature is only available on Model 74055. The active port is shown in the upper right corner of the display; “A” indicates the axial port (adjacent to the power switch) while “L” indicates the lateral port (opposite the input port).
- FILTE-** Move the 74010 Filter Wheel accessory to the specified position. Press FILTER, then type the number 1, 2, 3, 4, 5, or 6 indicating the desired position, then finally press ENTER to execute the move.
- SHIFT LABEL** This key (located directly above FILTER) provides access to the filter label. These labels are eight character alpha-numeric fields, used for information only. You can use the keypad to type any number, then press ENTER to accept that new value. The Hand Controller display only shows the first four digits of the label, and you can only enter numbers.
- LOCAL** Use this key to toggle between “local” communications through the Hand Controller and “remote” communication with a computer. Please note that if you press LOCAL to activate the Hand Controller (the display is lit and information fills the screen), you also lock out PC communications. You cannot command the MS260i™ via either RS-232, IEEE-488 or USB while the Hand Controller is active.
- SHIFT UNITS** To change the wavelength units press this key, followed by either 1, 2, or 3 and then press ENTER. Those three number keys are labeled with the corresponding units: 1 is nanometers (nm), 2 is micrometers ( $\mu\text{m}$ ), and 3 is wave numbers ( $\text{cm}^{-1}$ ) shown as “v”.
- 1** Pressing the 1 key when not in a function parameter edit mode brings up the “Slit1um>” prompt to edit/enter the slit 1 size, (the slit at the input port), the integer parameter is the open size specified in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6.
- SHIFT 1-** Pressing the SHIFT-1 key when not in a function parameter edit mode brings up the “Bandpass>” prompt. The floating point parameter is the wavelength resolution to automatically adjust slits for. If Bandpass is not set at zero the 74100 will be adjust all slits to maintain the Bandpass resolution at wavelength moves and grating changes. Setting Bandpass to 0 turns off this automatic adjustment. Slit range is 6-3000 microns, in 6 micron resolution steps. For Bandpass adjustments requiring slit openings larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6. Note; for Bandpass slit size calculations to function properly, Slit1 must be at the input port.
- 2** Pressing the 2 key when not in a function parameter edit mode brings up the “Slit2um>” prompt to edit/enter the slit 2 size, the integer parameter is the open size specified in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6.
- 3** Pressing the 3 key when not in a function parameter edit mode brings up the “Slit3um>” prompt to edit/enter the slit 3 size, the integer parameter is the open size specified in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6.

## IV.2 GETTING STARTED WITH IEEE-488

IEEE-488 communication (also called GPIB) offers the greatest versatility for integrating your MS260i™ into a larger system since you may connect up to 30 different devices to the same controller card. However, this versatility brings with it some complication in setting up each device.

Newport designs all of its IEEE-488 products to be completely compatible with National Instruments' GPIB controller cards. Likewise, most of our IEEE-488 software requires National Instruments' device drivers to run on your PC. We strongly recommend National Instruments' controller cards. Since there are differences between IEEE-488 implementations among the various manufactures, we cannot support them all. Therefore, Newport will only provide communication and software support for IEEE-488 systems that use National Instruments controller cards.

### Physical Connections

First, install your National Instruments control card and software in your PC computer. Run their diagnostic tools to make sure the card is working properly.

Next, be sure the MS260i™ power switch is off, then attach your IEEE-488 cable to the appropriate connector on the back panel. Attach the other end to the National Instruments' control card and secure the screws on both sides. Now turn on the MS260i™. In order to limit possible address conflicts, we recommend that you turn off all other IEEE-488 instruments while checking your MS260i™ for the first time. Note: the MS260i™ default GPIB primary address is 4.

### Verify Communications

You can check your IEEE-488 communications by running the application Cornerstone Utility Program . See software installation ( in section II. UNPACKING AND SETTING UP)

When this application runs, a dialog box will appear asking for the GPIB board index and Cornerstone™ 260 address. The default values, 0 and 4 respectively, should work for most systems. If you have multiple controller cards or if you have changed the Cornerstone™ 260 address fill in the correct values. Click the OK button to establish communication and open the main application window.

If you get an error at this point you may have a connection problem. Check all the cables (including the line cord), make sure the Cornerstone™ 260 is turned on, and make sure the Hand Controller is off. PLEASE NOTE: if the Hand Controller is active (the display is lit and information fills the screen), you cannot command the Cornerstone™ 260 via either RS-232 or IEEE-488. Press the LOCAL key on the Hand Controller to deactivate it, allowing PC communications. PLEASE ALSO NOTE: Cornerstone™ 260 selects either RS-232 or IEEE-488 communications based upon the first command received after power up. When you send an IEEE-488 command, or initialize your IEEE-488 software, the RS-232 link will be ignored until Cornerstone™ 260 power is turned off and on again.

### Low Level Communications

It may be desirable to send commands directly to the Cornerstone™ 260 through the National Instruments Interactive Control or through another simple interface. To establish low level communications with Cornerstone™ 260 you will have to first open the device. Using the National Instruments' drivers the correct command is: **ibdev 0 9 0 13 0 0x140a**

The timeout code "13" allows 30 seconds to read responses back from Cornerstone™ 260. In most cases 300 ms is sufficient (timeout code 10), however, the longer period guarantees success on all commands. The end of string code "0x140a" terminates a read when the linefeed character is

received from the instrument. All Cornerstone™ 260 responses end with carriage return and linefeed. Likewise, all messages sent to Cornerstone™ 260 must end with the linefeed character.

Again, refer to the detailed discussion of section V. FUNCTIONS AND CONTROLS for further information on direct low level communications, message syntax, and parameters.

### IV.3 GETTING STARTED WITH RS-232 COMMUNICATIONS

RS-232 connections are a simple means of communicating between your PC and your instruments. However, you are limited in that each instrument must have its own dedicated communication port at the computer. Most PC's come standard with COM1 and COM2, although COM1 is occasionally used by a serial mouse. Additional communication port hardware can be added and the corresponding drivers must be added to your Windows™ software. Please review your computer setup to determine what physical connections are available to you. Again, due to the variety of hardware and software options that are available Newport cannot provide support for PC configuration.

#### Physical Connections

First, make sure that the COM port you would like to use for MS260i™ is recognized by Windows™ and is working properly. Usually a simple check can be done with a serial mouse.

Next, be sure the MS260i™ power switch is off, then attach your RS-232 cable to the appropriate connector on the back panel. Attach the other end to the COM port connection on your PC. Note that RS-232 cables come in two types: modem and null-modem. The modem cable is a pin-to-pin cable. That is, each pin at one end connects directly to the corresponding pin at the other end: 1 to 1, 2 to 2, 3 to 3, etc. The modem cable is the correct cable for communicating with MS260i™.

Now turn on the MS260i™. You are ready to communicate with the instrument. The hardware communication parameters, including baud rate, are fixed inside MS260i™, so setup your COM port as follows:

Baud Rate:	9600
Data Bits:	8
Parity:	NONE
Stop Bits:	1
Flow Control:	NONE

#### Verify Communications

Check your RS-232 communications by running the Cornerstone Utility application. See software installation (in section II. UNPACKING AND SETTING UP).

When this application runs, a dialog box will appear asking for the COM port number. Fill in the correct value, then click the OK button to establish communication and open the main application window.

If you get an error at this point you may have a connection problem. Check all the cables (including the line cord), make sure the Cornerstone™ is turned on, and make sure the Hand Controller is off. PLEASE NOTE: if the Hand Controller is active (the display is lit and information fills the screen), you cannot command the Cornerstone™ via either RS-232 or IEEE-488. Press the LOCAL key on the Hand Controller to deactivate it, allowing PC communications. PLEASE ALSO NOTE: Cornerstone™

selects either RS-232 or IEEE-488 communications based upon the first command received after power up. When you send an RS-232 command, or initialize your RS-232 software, the IEEE-488 link will be ignored until Cornerstone™ power is turned off and on again.

### **Low Level Communications**

In some cases, it may be desirable to send commands directly to Cornerstone™ through a terminal emulation program, such as the HyperTerminal™ provided with Windows™. Once your terminal program is open and running you will have to first select a COM port and configure it according to the parameters previously listed (9600, 8, N, 1, None). Additionally, you should set the terminal to "Send line ends with line feeds". There is no reason to turn on the local echo since Cornerstone™ will automatically echo each character as it is typed. All Cornerstone™ responses end with carriage return and linefeed. Likewise, all messages sent to Cornerstone™ must end with the linefeed character.

Again, refer to the detailed discussion of section V. FUNCTIONS AND CONTROLS for further information on direct low level communications, message syntax, and parameters.

---

## V. FUNCTIONS AND CONTROLS

---

### V.1 COMMAND AND QUERY SYNTAX

For the purposes of this discussion we will divide communications with the instrument into two parts. First, the messages that you send to MS260i™ will be called “Statements”. Second, the messages that you receive from MS260i™ will be called “Responses”. When communicating with MS260i™ you will use two types of Statements: “Commands” and “Queries”. A Command causes some physical action or sets an internal parameter. Queries ask some question of the instrument and automatically return a Response. Fundamentally, the syntax is the same for all messages.

#### Termination

Both Statements and Responses end with a carriage return (ASCII character code 13 decimal) and linefeed (ASCII character code 10 decimal). This manual will show the termination as [cr][lf]. Please note, you may find it more convenient to end Statements with only the linefeed [lf]; this is also acceptable.

#### Capitalization

Statements may be sent in upper case, in lower case, or in any combination of the two. Responses will always be returned in upper case. For clarity throughout this manual, Statements will be shown in upper case (ie. STATEMENT) and Responses will be shown in upper case italics (ie. *RESPONSE*).

#### Parameters

Some Statements will require a parameter. In this manual numeric parameters will be shown as “XX”; each X will fill a numeric position and decimals are used to indicate where floating point number should be expected. Numeric notation is not strict and integers can be substituted for floating point numbers as desired. There are some specific cases where alpha-numeric values or exponential notation may be required; these special Statements will be detailed in the sections that follow.

Most normal communication with MS260i™, especially communication handled inside closed application software, is done in Standard Mode. This mode follows all of the rules on the previous page plus there are some additional features for error handling.

MS260i™ stores errors internally in a status byte and an error byte, where each bit has a specific meaning. Details of the status and error bytes can be found in this manual under section V-4 ERROR CODES. For now, we will simply cover the syntax necessary to send Statements and receive Responses in this communication mode.

In all cases, Statements must be terminated properly with [lf] for MS260i™ to take any action. Statements that are not understood or parameters that are unacceptable will generate an error code in the status byte, ie. status byte ≠ 0. Moreover, the status byte must be explicitly read, it is not returned after general Statements. Correct Statements are handled as follows:

#### Commands

Under RS-232, Statements are echoed back to the PC. Thus, after sending a Command you should read back this echo. There is no further response to a Command.

For example:            Send: GOWAVE 500[lf]            Receive: *GOWAVE 500[lf]*

Under IEEE-488, there is no echo. There is no automatic response to a Command.

For example:            Send: GOWAVE 500[lf]            Receive: {none}

**Queries**

Under RS-232, Statements are echoed back to the PC. Additionally, there will be an explicit response to that query. Thus, after sending a Query you should read back the echo and the answer together.

For example:      Send : WAVE?[lf]      Receive: *WAVE?[lf]500[cr][lf]*

Under IEEE-488, there is no echo. Thus, you can read back the Response by itself.

For example:      Send : WAVE?[lf]      Receive: *500[cr][lf]*

**V.2 HANDSHAKING MODE**

This communications mode is most useful for methods development, troubleshooting, or in the early stages of programming. Use the command: "HANDSHAKE 1[lf]" to put MS260i™ into Handshake Mode. This mode follows the command syntax listed at the start of section V-1 in this manual, but in addition every Statement is acknowledged. The "handshake" that you receive is the status byte.

MS260i™ stores errors internally in a status byte and an error byte, where each bit has a specific meaning. Details of the status and error bytes can be found in this manual under section V-4 ERROR CODES. For now, we will simply cover the syntax necessary to send Statements and receive Responses.

In all cases, Statements must be terminated properly with [lf] for MS260i™ to take any action. Statements that are not understood or parameters that are unacceptable will generate an error code in the status byte, ie. status byte  $\neq$  0. This status byte will be reported immediately if it contains an error, otherwise the status byte "00" will be returned upon completion of the Statement. Correct Statements are handled as follows:

**Commands**

Under RS-232, Statements are echoed back to the PC. Thus, after sending a Command you should read back this echo. Additionally, the status byte is reported when the action is complete.

For example:      Send: GOWAVE 500[lf]      Response: *GOWAVE 500[lf]00[cr][lf]*

Under IEEE-488, there is no echo. The only response will be the status byte.

For example:      Send: GOWAVE 500[lf]      Response: *00[cr][lf]*

**Queries**

Under RS-232, Statements are echoed back to the PC. Additionally, there will be an explicit response to that Query, and the status byte. Thus, after sending a Query you should read back a 3 part answer.

For example:      Send : WAVE?[lf]      Response: *WAVE?[lf]500[cr][lf]00[cr][lf]*

Under IEEE-488, there is no echo. Thus, you will read back the Response and the status byte.

For example:      Send : WAVE?[lf]      Response: *500[cr][lf]00[cr][lf]*



**V.3 BRIEF COMMAND REFERENCE ORGANIZED ALPHABETICALLY**

<b>Command</b>	<b>Query</b>	<b>Hand Ctrl Key</b>
ABORT		[ABORT]
ADDRESS XX	ADDRESS?	
CALIBRATE XXX.XXX		[CALIB]
	ERROR?	
FILTER X	FILTER?	[FILTER]
FILTER1LABEL TTTTTTTT	FILTER1LABEL?	[SHIFT LABEL]
GOWAVE XXX.XXX	WAVE?	[GO WAVE]
GRAT X	GRAT?	[GRAT]
GRAT1FACTOR XXX.XXX	GRAT1FACTOR?	
GRAT1LABEL TTTTTTTT	GRAT1LABEL?	[SHIFT LABEL]
GRAT1LINES XXXX	GRAT1LINES?	[LINES]
GRAT1OFFSET XXX.XXX	GRAT1OFFSET?	
HANDSHAKE X	HANDSHAKE?	
INFO?		[LOCAL]
OUTPORT T	OUTPORT?	[PORT]
SHUTTER T	SHUTTER?	[SHUTTER]
	STB?	
STEP XXX	STEP?	[STEP UP]
UNITS TT		[SHIFT UNITS]
SLIT1MICRONS XXXX	SLIT1MICRONS?	[1]
SLIT2MICRONS XXXX	SLIT2MICRONS?	[2]
SLIT3MICRONS XXXX	SLIT3MICRONS?	[3]
BANDPASS XXX.XX	BANDPASS?	[SHIFT 1]

#### V.4 DETAILED COMMAND REFERENCE ORGANIZED BY FUNCTION

Recall from section V-1 of this manual that messages that you send to MS260i™ will be called “Statements” and messages that you receive from MS260i™ will be called “Responses”. There are two types of Statements: “Commands” and “Queries”. A Command causes some physical action or sets an internal parameter. Queries ask some question of the instrument and automatically return a Response. Also recall that all Statements must end in [lf], and [cr][lf] is acceptable. Likewise, all Responses will end with [cr][lf]. Refer to section V-1 for detailed general syntax instructions. This section describes the exact syntax, acceptable parameters, and functionality for every MS260i™ Command. For each corresponding Query, in braces {}, we also list the type of Response you should expect. Where appropriate the corresponding Hand Controller key, in brackets [ ], is also listed. Finally, there are some warnings as to possible errors that could occur if a command is not used properly.

##### MS260i™ Motion Commands:

UNITS TT {UNITS?} [UNITS]  
 Specifies the operational unit designated by the mnemonic parameter “TT”. Possible parameters are: “NM” for nanometers “UM” for micrometers “WN” for wavenumbers (cm<sup>-1</sup>)

By default MS260i™ uses nanometers. Parameters for all motion Statements are handled in the current units. Therefore, if you change from nanometers to micrometers GOWAVE parameters must be entered in micrometers. Likewise, WAVE? answers will be returned in micrometers. The response to UNITS? is always the two character mnemonic for the current units.

GOWAVE XXX.XXX {WAVE?} [GO WAVE]  
 Moves the wavelength drive at maximum speed to the step position closest to the specified wavelength parameter. This parameter must be in the current units and may be either a floating point number or an integer, ie. “GOWAVE 500.00[lf]” and “GOWAVE 500[lf]” will be understood the same way. For each grating there is a maximum possible wavelength (approximately 1600 nm for a 1200 l/mm grating, thus 800 nm for a 600 l/mm, etc.) This command will generate an error if the parameter is larger than the maximum wavelength. The response to WAVE? will be a floating point number indicating the exact wavelength position. Current units are assumed and so no units are passed with the floating point number.

CALIBRATE XXX.XXX [CALIB]  
 This command is used to make minor corrections to the wavelength calibration of MS260i™. It is used when installing new gratings, or at any point when the wavelength calibration is disrupted. Execute this command to define the current position as the wavelength specified in the numeric parameter XXX.XXX. This calculates a new OFFSET for the grating and saves it accordingly. Please Note: The calibration is corrected for the entire range of the current grating and current output port; on a 74055 system, a separate calibration may be necessary if the output port is changed. A complete calibration procedure is outlined in Section VII-4 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION.

ABORT [ABORT]  
 Simply stops any wavelength motion immediately, ie. GOSCAN, GOWAVE, STEP, etc.

STEP XXXX {STEP?} [STEP UP] / [STEP DOWN]  
 Move the wavelength drive by the integer number of steps indicated in XXXX. The parameter can be positive or negative, however, one step is the smallest increment MS260i™ can move and so this parameter must be an integer. The response to STEP? is the current step position as measured from

the physical home sensors.

### MS260i™ Grating Commands:

**GRAT X** {GRAT?} [GRAT]  
 Move the wavelength drive to the grating specified in the integer parameter. The parameter should be either "1", "2" or "3". Please note: whenever you change gratings, the wavelength drive will sweep past zero order on grating #2. In order to prevent saturation of a detection system, it is wise to close the shutter temporarily while changing gratings. When changing to grating #1 the instrument stops at the maximum wavelength for that grating. When changing to grating #3, the system stops at the zero order position for that grating. When changing to grating #2, the system will stop at the zero order position or the maximum wavelength, whichever is closer. Therefore, this command is usually followed by a GOWAVE command. The response to GRAT? contains three parameters separated by commas without spaces. The first parameters is the grating number, then the lines/mm, then the label for that grating. For example: "1,1200,BLUE[cr][lf]".

**GRAT1LABEL TTTTTT** {GRAT1LABEL?} [SHIFT LABEL]  
 Sets the user defined "label" for the grating #1 to the < 9 character parameter "TTTTTTT". The parameter can letters or numbers. This label is for informational purposes only. Newport uses these labels to indicate blaze wavelength for those gratings that are installed in MS260i™ at the factory. Note: that the Hand Controller only displays four characters, however all eight may be entered. Also, the Hand Controller only changes the current grating parameters.

**GRAT2LABEL TTTTTT** {GRAT2LABEL?} [SHIFT LABEL]  
 Sets the user defined "label" for the grating #2 to the < 9 character string parameter "TTTTTTTT". The parameter can letters or numbers. This label is for informational purposes only. Newport uses these labels to indicate blaze wavelength for those gratings that are installed in MS260i™ at the factory. Note: that the Hand Controller only displays four characters, however all eight may be entered. Also, the Hand Controller only changes the current grating parameters.

**GRAT3LABEL TTTTTT** {GRAT3LABEL?} [SHIFT LABEL]  
 Sets the user defined "label" for the grating #3 to the < 9 character string parameter "TTTTTTTT". The parameter can letters or numbers. This label is for informational purposes only. Newport uses these labels to indicate blaze wavelength for those gratings that are installed in MS260i™ at the factory. Note: that the Hand Controller only displays four characters, however all eight may be entered. Also, the Hand Controller only changes the current grating parameters.

**GRAT1LINES XXXX** {GRAT1LINES?} [LINES]  
 Sets the lines per millimeter of grating #1 to the integer parameter specified as XXXX. Warning: this value is used to calculate wavelength motion. Line/mm are usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. Note, the Hand Controller only changes the current grating parameters.

**GRAT2LINES XXXX** {GRAT2LINES?} [LINES]  
 Sets the lines per millimeter of grating #2 to the integer parameter specified as XXXX. Warning: this value is used to calculate wavelength motion. Line/mm are usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. Note, the Hand Controller only changes the current grating parameters.

**GRAT3LINES XXXX** {GRAT3LINES?} [LINES]

Sets the lines per millimeter of grating #3 to the integer parameter specified as XXXX. Warning: this value is used to calculate wavelength motion. Line/mm are usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. Note, the Hand Controller only changes the current grating parameters.

GRAT1FACTOR XXX.XXX {GRAT1FACTOR?}

Sets the calibration factor of grating #1 to the numeric parameter specified as XXX.XXX. Warning: this value is used to calculate wavelength motion. Factor is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport.

GRAT2FACTOR XXX.XXX {GRAT2FACTOR?}

Sets the calibration factor of grating #2 to the numeric parameter specified as XXX.XXX. Warning: this value is used to calculate wavelength motion. Factor is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport.

GRAT3FACTOR XXX.XXX {GRAT3FACTOR?}

Sets the calibration factor of grating #3 to the numeric parameter specified as XXX.XXX. Warning: this value is used to calculate wavelength motion. Factor is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport.

GRAT1OFFSET XXX.XXX {GRAT1OFFSET?}

Sets the calibration offset for grating #1 as the floating point parameter specified in XXX.XXX. Warning: this value is used to calculate wavelength motion. Offset is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. This value is automatically changed via the CALIBRATE command. A calibration procedure is outlined in Section VII-4 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION.

GRAT2OFFSET XXX.XXX {GRAT2OFFSET?}

Sets the calibration offset for grating #2 as the floating point parameter specified in XXX.XXX. Warning: this value is used to calculate wavelength motion. Offset is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. This value is automatically changed via the CALIBRATE command. A calibration procedure is outlined in Section VII-4 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION.

GRAT3OFFSET XXX.XXX {GRAT3OFFSET?}

Sets the calibration offset for grating #3 as the floating point parameter specified in XXX.XXX. Warning: this value is used to calculate wavelength motion. Offset is usually changed only when installing a new grating, and in that case the parameter should match the value provided by Newport. This value is automatically changed via the CALIBRATE command. A calibration procedure is outlined in Section VII-4 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION.

### MS260i™ Accessory Commands:

SHUTTER T {SHUTTER?} [SHUTTER]

Closes or opens the shutter depending upon the single letter mnemonic parameter "T". Closing the shutter requires "C", whereas opening it requires "O". The response to SHUTTER? is simply the one character parameter describing the current state. When using the Hand Controller key [SHUTTER] no parameter is necessary. Pressing this key toggles the shutter out of the current state into the other state. Look to the display to check shutter position.

OUTPORT T                      {OUTPORT?}                      [OUTPORT]

74055 has two output ports; this command moves the motorized flip mirror to activate the output port indicated by the single letter mnemonic parameter (T). Use "1" to select the axial port (default and adjacent to the power switch) and "2" to select the lateral port. The response to PORT? Is simply the one character parameter describing the current state. When using the Hand Controller key [PORT] no parameter is necessary. Pressing this key toggles the flip mirror out of the current state into the opposite position. Look to the top right corner of the display to check the mirror position.

FILTER X                      {FILTER?}                      [FILTER]

Moves the filter wheel to the position specified in the single integer parameter "X". Acceptable values range from 1 through 6. The response to FILTER? is the integer value describing the current filter position; if there is no filter wheel connected, or if the wheel is out of position, "0" is returned and an error is set in the status byte. Sending this command while the filter wheel is out of position will cause the wheel to search for #1 and then move to the requested position. This search occurs at power on too.

FILTER#LABEL TTTTTTTT {FILTER#LABEL?}                      [SHIFT LABEL]

Sets the label for the filter position specified by # to the < 9 character string parameter "TTTTTTTT". Valid values of # are 1-6, using any other value will result in a status byte error indicating a bad command, not a bad parameter. The value of "TTTTTTTT" can be letters or numbers. This label is for informational purposes only.

SLIT1MICRONS XXXX                      {SLIT1MICRONS?}                      [1]

Commands the MS260 to open slit 1, (the slit at the input port), to the size specified by the integer parameter in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6. If the query responds with 0, there is a problem with the slit.

SLIT2MICRONS XXXX                      {SLIT2MICRONS?}                      [2]

Commands the MS260 to open slit 2 to the size specified by the integer parameter in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6. If the query responds with 0, there is a problem with the slit.

SLIT3MICRONS XXXX                      {SLIT3MICRONS?}                      [3]

Commands the MS260 to open slit 3 to the size specified by the integer parameter in microns. Slit range is 6-3000 microns, in 6 micron resolution steps; for values larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6. If the query responds with 0, there is a problem with the slit.

BANDPASS                      XXX.XX                      {BANDPASS?}                      [SHIFT 1]

The MS260 uses the Bandpass parameter to automatically adjust all the slits. The floating point parameter is the wavelength resolution to automatically adjust slits for. If Bandpass is not set at zero the 74000 will be adjust all slits to maintain the Bandpass resolution at wavelength moves and grating changes. Setting Bandpass to 0 turns off this automatic adjustment. Slit range is 6-3000 microns, in 6 micron resolution steps. For Bandpass adjustments requiring slit openings larger than 3000 the slit goes to 3000, for values 0-6 the slit goes to 6. Note; for Bandpass slit size calculations to function properly, Slit1 must be at the input port.

**MS260i™ System Commands:**

{INFO?} [LOCAL]

Query for basic instrument information. Generates the same response as the GPIB specific \*IDN? That response is "Newport, Model 74050 MS260i™,SNXXX,VYY.YY.YY" where XXX is the unit's serial number and YY.YY.YY is the version number for internal firmware. This information is displayed for 3 seconds when [LOCAL] is pressed on the Hand Controller.

HANDSHAKE X {HANDSHAKE?}

This command turns Handshake Mode on and off. Setting the parameter "X" to "1" causes MS260i™ to acknowledge command completion according to the description given in Section V-1-B. HANDSHAKE MODE. Setting the parameter to "0" returns MS260i™ to normal operation according to the description given in Section V-1-A STANDARD MODE. The response to the query HANDSHAKE? is the single integer describing the current mode, "0" or "1".

{STB?}

Use this query to check for MS260i™ error conditions. The response will be "32" when an error occurred or "00" when no error is present. Reading this code clears the status byte. Further information can be obtained using the ERROR? query below. A detailed discussion of error handling is given in section V-4.{ERROR?}

Use this query to obtain further information about an error state reported in the Status byte. The response will be an integer from 0 through 9. The definition of each error code and a detailed discussion of error handling is given in section V-4.

ADDRESS XX {ADDRESS?}

Use this command to change the IEEE-488 primary address for MS260i™ if the default address (9) conflicts with another instrument in your laboratory. You should turn off the other instrument, establish communication with MS260i™, and send this command with the integer parameter from 1 to 30. If successful, the address change takes place immediately and any further communication using the old address will cause a IEEE-488 error. If unsuccessful an error will be reported in the Status byte.



## VI. MS260i™ UTILITY PROGRAM

MS260 software is loaded onto your PC per the directions in Section II UNPACKING AND SETTING UP. The Utility Program comes with three versions: IEEE-488, RS-232 and USB communications. This program provides easy access to most of Cornerstone™'s functions. It is especially useful for adding new gratings to your instrument; follow the instructions in Section VII-4 ADDITIONAL GRATINGS - WAVELENGTH CALIBRATION to add new gratings.

### Main Window (Refer to Figure 1, Page 25)

The "Present Wavelength" indicates the wavelength position of the monochromator. All the wavelength units indicators on the Utility program's main screen are determined by the setting the Units Dialog Box.

The VCR type buttons under the "Present Wavelength" field can be used to position the monochromator by steps; reverse slew, back one step, stop, forward step, and fast forward.

The numeric control beneath the step motion buttons is used to position the monochromator by wavelength; enter the desired wavelength position and click on the "GoTo" button.

The Shutter status indicator is "Shutter Open"/"Shutter Closed."

The "Grating" indicator shows the active grating and its lines/mm.

The "Filter" indicator shows the present filter wheel position.

The "Scan Parameters" section enables the user to edit scan range, interval and time for the "Go Scan" button.

"Start Wavelength"	scan to begin at this wavelength
"End Wavelength"	scan to end at this wavelength
"Interval Wavelength"	scan to pause at interval wavelength points
"Pause"	seconds to wait at each scan point

Clicking "Go Scan" will run a monochromator scan according to the "Scan Parameters"

The buttons on the right side of the Utility Program's main window will bring up dialog boxes for different functions:

"Units"	Used to select the monochromator wavelength units, nanometers, micrometers or wave numbers.
"Calibrate"	Bring up a dialog to allow the user to calibrate the monochromator by entering the actual wavelength position.
"Setup"	Allows the editing/display of the monochromator's grating parameters, should never be needed unless installing new gratings.
"Shutter"	Allows control of the shutter.
"Grating"	Allows selection of a specific grating, or enables automatic grating selection for specific wavelength regions.
"Filter"	Allows selection of a specific filter wheel position, or enables automatic filter position for specific wavelength regions.



- “Setup Comm” Used to select IEEE-488 or RS-232 communication parameters, i.e. board address or port number. Not needed for USB (74125).
- “Exit” Exits the utility program and saves monochromator and communication parameters.

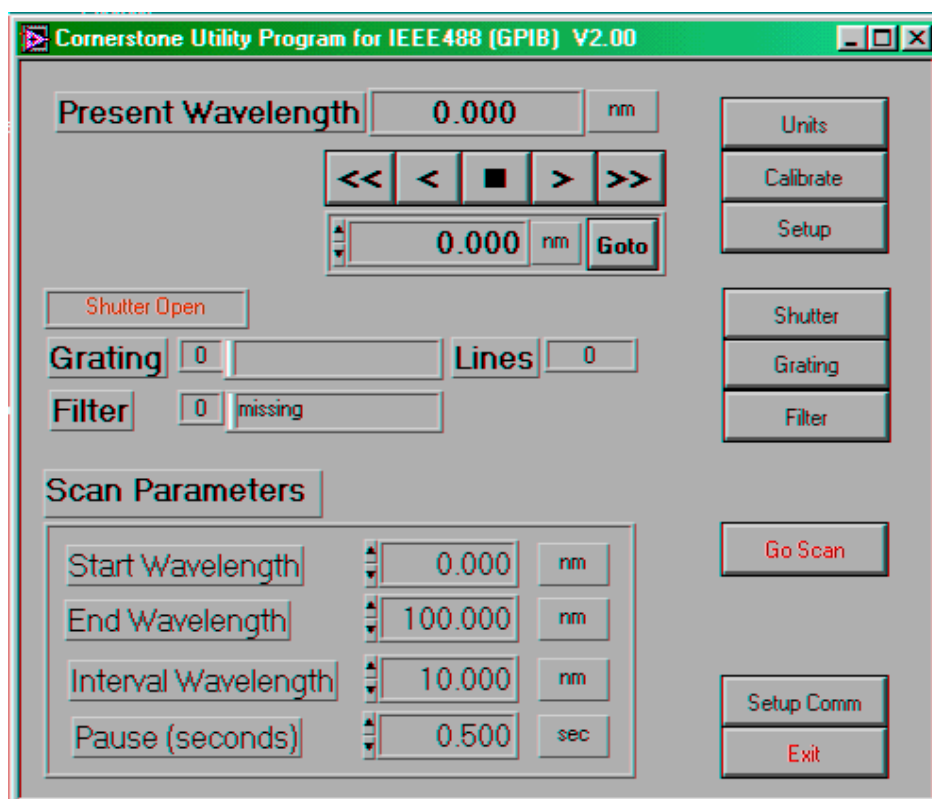


Figure 1: MS260 Utility Program, Main Window

## VI.1 SOURCE CODE

This program is a closed executable file written in National Instrument's LabView™ and is based on our collection of 74085/74086™ VI's. The Utility Program diagram may serve as an example or starting point for further development in LabView™. If you are a LabView™ programmer and own a full LabView™ development environment, you can load the VI's from your hard drive. There are three versions of the VI's, one each for RS-232, IEEE-488 and USB. See in Section II Table I for the location of the VIs on the CD.

## VI.2 IEEE-488 VERSION

This version of the Utility Program requires that you have a National Instrument's IEEE-488 Controller Card and associated software drivers installed in your PC. It is important to configure your PC and establish communications before proceeding with any application software. Please follow the procedure in Section IV-2 GETTING STARTED WITH IEEE-488 COMMUNICATION to check your communications before running this Utility Software.

The program will run and a dialog box will appear asking for the GPIB board index and Cornerstone™ address. The default values, 0 and 4 respectively, should work for most systems. If you had to change these values to run the diagnostic software in Section IV-2, fill in the correct values here again. Click the OK button to establish communication and open the main application window.

### **VI.3 RS-232 VERSION**

Again in this case, it is important to configure your PC and establish communications before proceeding with any application software. Please follow the procedure in Section IV-3 GETTING STARTED WITH RS-232 COMMUNICATION to check your communications before running this Utility Software.

The program will run and a dialog box will appear asking for the COM port number. Fill in the correct value, then click the OK button to establish communication and open the main application window.

### **VI.4 USB VERSION 74085/74086**

A USB2 port must be available on the PC. The ORIEL\_CS\_USB driver must be installed before the 74125 can be controlled via a PC. Please see Section II on Software Installation.

---

## VII. OPTIONAL ACCESSORIES

---

### VII.1 MICROMETER DRIVEN SLIT

The 74001 slit is adjustable from approximately 4  $\mu\text{m}$  to 3 mm using the micrometer drive on top. You should apply a 10x multiplier to the micrometer markings to readout the slit opening. Thus, each mark on the micrometer barrel corresponds to 10  $\mu\text{m}$  of slit width. The height is also continuously adjustable from 2 mm to 15 mm tall. Please note that optimum spectral resolution for any monochromator is obtained with short, narrow slits. However, the 74001 is designed primarily for versatility and convenience in changing resolution and throughput; fixed slits should be used for the utmost accuracy and repeatability, especially at high resolution.

Slit Width:	4 $\mu\text{m}$ to 3 mm
Slit Height:	2 mm to 15 mm
Precision:	+/- 10 $\mu\text{m}$
Accuracy:	+/- 10 $\mu\text{m}$ to 250 $\mu\text{m}$
	+/- 5% from 250 $\mu\text{m}$ to 3 mm

### VII.2 FILTER WHEEL

Newport's model 74010 Motorized Filter Wheel is matched for use with MS260i™. The drive electronics and command set for this wheel are included with every MS260i™ and the wheel mounts directly to the monochromator via Newport's standard 1.5" Series Flange. The 74010 is a six position enclosed wheel that holds 1 inch (25 mm) diameter filters, with a clear aperture of 0.8 inches (20 mm). Each filter is mounted in its own holder and the holders are individually removable. Moreover, when you change a filter, there is no need to break your optical path since the filter holders can be exchanged from the top of the wheel.

Please note: None of the MS260i™ accessories are designed to be connected or disconnected with power on. Always turn off the instrument using the power switch on the back panel before you connect or remove the filter wheel.

6 position Motorized Filter Wheel	Model 74010
1" Filter Holder (for use in wheel)	Model 74011
Rod Mount for 74011 Holders	Model 74012
1.5" Series Mount for 74011	Model 74013

### VII.3 ADDITIONAL GRATINGS- WAVELENGTH CALIBRATION

Most MS260i™ Spectrographs are shipped with three gratings installed. However, these gratings are individually replaceable. That is, a single grating can be removed and a new one mounted in its place at your laboratory. There is no need to purchase more than one grating or a larger grating turret, and there is no need to send the instrument back to Newport. We offer a wide variety of gratings to suit most spectral ranges and resolution requirements. Please call Newport with your specific needs.

#### Installation

Each grating comes mounted and aligned on a completely adjustable grating mount. For most applications there is no need to adjust this grating holder. Simply follow the steps below to mount your new grating:

1. Using the Hand Controller, or directly from the PC, close the shutter and command MS260i™ to go to grating #1. Then send the instrument to zero order. (GOWAVE 0).

2. Turn off the power switch and unplug the instrument.
3. Remove the cover screws and lift it off. (WARNING: the optics inside this instrument are coating to protect and enhance their reflectivity. This coating is extremely soft and easily damaged. Do not touch any of the optics. Do not attempt to clean any of the optics, except with a gentle stream of filtered dry nitrogen. Since accidents do happen, we recommend wearing powder free, latex gloves when working inside the optical compartment of any instrument.)
4. There are three long screws, one on each side of the gratings. Grab the knurled top with you fingers, loosen and remove each screw.
5. The grating that is facing the main mirrors is grating #1, while the grating that faces the electronics compartment is grating #2. Only change one grating at a time. Decide which one you would like to remove, grab the sides of that grating holder with one hand, and firmly pull the grating up and out of the instrument. Be careful not to remove the other gratings; you may want to hold it in place using one finger from your other hand.
6. Uncover your new grating, and use that cover to protect the grating you have just removed.
7. Note that inside the instrument, extending up from the top of the brass gear where the grating sits, is a small aluminum pin. This pin will locate your new grating.
8. Slide the new grating into the instrument. The pin mentioned in step #7 should come up through the clearance hole in the bottom of the grating mount. Make sure that the grating mount sits flat on the grating base.
9. If you are replacing more than one grating, return to steps #5 - #8 for the next grating.
10. Now replace the long knurled screws that hold the gratings in place. It is best to alternate between the screws, not to tighten one completely and then the next. The screws should line up parallel with the grating face. These screws do not require excessive pressure to hold the gratings in place, use only
11. firm finger pressure to tighten them.
12. Replace the cover and secure the screws. Again, start all of the screws and then tighten each one a little at a time- do not completely tighten one and then start the next.
13. Plug in the power cable and turn on the power switch.
14. Run the MS260i™ Utility Program of your choice, either IEEE-488 or RS-232. If you have not run this software before, refer to the instructions in Section VI. CORNERSTONE UTILITY PROGRAM.
15. Once the software is running and you see the main window, press the Grating Parameters button in the center of the control panel.
16. Now enter you new grating parameters, as provided by Newport, in the correct column.
17. Press SEND at the bottom of the screen. Your grating is now installed and updated.

**Grating Factor Calibration**

1. Select a radiation source that has at least two narrow spectral lines in the wavelength region where you will be working.
2. The radiation from your line source should be focused on the entrance slit to fill the monochromator's f/3.7.
3. The focused beam should be parallel to Cornerstone™'s optical axis.
4. Command Cornerstone™ to move to the grating you would like to calibrate and then to a wavelength of your line source. Now use the STEP feature to move Cornerstone™ toward the signal peak for your source.
5. Notate the spectral line you first commanded to move to and the resultant wavelength position where you located the actual peak. See the tables below.
6. Command Cornerstone™ to move to a second wavelength of your line source. Now use the STEP feature to move Cornerstone™ toward the signal peak for your source.
7. Notate the spectral line you first commanded to move to and the resultant wavelength position where you located the actual peak. See the example tables below.
8. Using Factor equation below to calculate these parameters for your new grating.
9. Repeat for all gratings, Using the Utility software you can enter the FACTOR values into the Cornerstone via the "Setup" screen. The GRATINGnFACTOR commands can be used, VIs are also supplied for these commands. They cannot be entered directly via the Hand Controller.

	Spectral Line	Observed Peak
Wavelength 1	L1	P1
Wavelength 2	L2	P2

$$\text{FACTOR} = |L1 - L2| / |P1 - P2|$$

**EXAMPLE:**

	Spectral Line	Observed Peak
Wavelength 1	546	577
Wavelength 2	365	372

$$\text{Factor} = |577 - 365| / |577 - 372| = 181 / 205 = 0.8829268$$

Blank tables to notate your peak locations

**GRATING1**

	Spectral Line	Observed Peak
Wavelength 1		
Wavelength 2		

GRATING 1 FACTOR =

## GRATING 2

	Spectral Line	Observed Peak
Wavelength 1		
Wavelength 2		

GRATING 2 FACTOR =

## GRATING 3

	Spectral Line	Observed Peak
Wavelength 1		
Wavelength 2		

GRATING 3 FACTOR =

## Wavelength Calibration

The Cornerstone comes calibrated from the factory you can use the Cornerstone CALIBRATE command to adjust offset, this can be done via PC or with the Hand Controller.

For the highest accuracies, it may be necessary under some circumstances to wavelength calibrate your MS260i™. We recommend using a low pressure spectral line lamp such as Newport's model 6035 Mercury Lamp for this purpose, however you can use another line source such as a laboratory laser. Set up for this task is quite important. Follow these guidelines:

1. Select a radiation source that has at least one narrow spectral line in the wavelength region where you will be working.
2. The radiation from your line source should be focused on the entrance slit to fill the monochromator's f/3.7.
3. The focused beam should be parallel to MS260i™'s optical axis.

Once you have set up the known source and your detection system, wavelength calibration is quite simple. First, command MS260i™ to move to the grating you would like to calibrate and then to the wavelength of your line source. Second, use the STEP feature to move MS260i™ toward the signal peak for your source. Finally, when you have placed the peak in the center of your CCD, use the CALIBRATE feature (CALIB on the Hand Controller) to enter the exact wavelength of your spectral line. MS260i™ will recalculate its motion to place that wavelength in the position you have specified, and it will correct all other wavelength positions as well. However, you may need to repeat this for the another grating, or if you add a new grating later.

This chapter suggests solutions to potential problems you may encounter when operating Cornerstone. This section will be periodically updated on our web site. If you have a question that is

not answered in either place, please call a Newport Applications Engineer for assistance.

#### **VII.4 MEMORY LOSS (LARGE WAVELENGTH CALIBRATION ERRORS)**

**Symptoms:** 1- there are large errors in wavelength positioning of all gratings (maybe 100nm).  
2- the instrument lists all gratings as 1200 l/mm.

**Problem:** The instrument memory has been cleared. This may be the result of a power failure or a spike on the mains. It may also result from electro-static discharge or other high field radiation in contact with, or very nearby, you instrument.

**Solution:** You will have to reload the instrument parameters. All of the necessary values are included on the Cornerstone Calibration Parameters datasheet that came with your instrument. First, run the Utility program that you installed from the Cornerstone CD (according to section II of this manual). You can enter the calibration parameters via the Utility programs "Setup" screen.

**VIII. DECLARATION OF CONFORMITY**

**Manufacturer's Name:** Newport Corporation

**Manufacturer's Address:** 150 Long Beach Boulevard  
Stratford, CT 06615  
USA

**Declares that the product**

**Product Name:** MS260i™ 1/4m Imaging Spectrograph

**Model Number:** 74050, 74055, 74085 and 74086

**Conforms to the following Product Specifications:**

**Safety:** EN 61010-1: 1993 + A2/ IEC 1010-1: 1990 + A1 + A2

**EMC:** EN 50081-1: 1992  
EN 55022: 1994/ EN 55011: 1993 Class B

EN 50082-1: 1992  
IEC 801-2: 1991/ IEC 1000-4-2: 1995  
ENV 50140:1993/ EC 1000-4-3:1995/ EN 61000-4-3: 1995  
IEC 801-4: 1988/ IEC 1000-4-4: 1995

**Complies with the following Directives:**

- the EMC Directive 89/336/EEC

- the Low Voltage Directive 73/23/EEC

**And, accordingly, carries the CE mark.**

Stratford, CT June, 1998



George Buzel

Director of Engineering



## IX. WARRANTY AND RETURNS

Newport warrants that all goods described in this manual (except consumables such as lamps, bulbs, filters, ellipses, etc.) shall be free from defects in material and workmanship. Such defects become apparent within the following period:

1. All products described here, except spare parts: one (1) year or 3000 hours of operation, whichever comes first, after delivery of the goods to the buyer.
2. Spare parts: ninety (90) days after delivery of goods to the buyer.

Newport's liability under this warranty is limited to the adjustment, repair and/or replacement of the defective part(s). During the above listed warranty period, Newport shall provide all materials to accomplish the repaired adjustment, repair or replacement. Newport shall provide the labor required during the above listed warranty period to adjust, repair and/or replace the defective goods at no cost to the buyer ONLY IF the defective goods are returned, freight prepaid, to a Newport designated facility. If goods are not returned to Newport, and the user chooses to have repairs made at their premises, Newport shall provide labor for field adjustment, repair and/or replacement at prevailing rates for field service, on a portal-to-portal basis.

Newport shall be relieved of all obligations and liability under this warranty of:

1. The user operates the device with any accessory, equipment or part not specifically approved or manufactured or specified by Newport unless buyer furnishes reasonable evidence that such installations were not the cause of the defect. This provision shall not apply to any accessory, equipment or part which does not affect the safe operation of the device.
2. The goods are not operated or maintained in accordance with Newport's instructions and specifications.
3. The goods have been repaired, altered or modified by other than authorized Newport personnel.
4. Buyer does not return the defective goods, freight prepaid, to a Newport facility within the applicable warranty period.

IT IS EXPRESSLY AGREED THAT THIS WARRANTY SHALL REPLACE ALL WARRANTIES OF FITNESS AND MERCHANTABILITY. BUYER HEREBY WAIVES ALL OTHER WARRANTIES, GUARANTEES, CONDITIONS OR LIABILITIES, EXPRESSED OR IMPLIED, ARISING BY LAW OR OTHERWISE, WHETHER OR NOT OCCASIONED BY NEWPORT'S NEGLIGENCE.

This warranty shall not be extended, altered or varied except by a written document signed by both parties. If any portion of this agreement is invalidated, the remainder of the agreement shall remain in full force and effect.

### CONSEQUENTIAL DAMAGES

Newport shall not be responsible for consequential damages resulting from misfunctions or malfunctions of the goods described in this manual. Newport's total responsibility is limited to repairing or replacing the malfunctioning or malfunctioning goods under the terms and conditions of the above described warranty.

### INSURANCE

Persons receiving goods for demonstrations, demo loan, temporary use or in any manner in which title is not transferred from Newport, shall assume full responsibility for any and all damage while in their care, custody and control. If damage occurs, unrelated to the proper and warranted use and performance of the goods, recipient of the goods accepts full responsibility for restoring the goods to their condition upon original delivery, and for assuming all costs and charges.

### RETURNS

Before returning equipment to Newport for repair, please call the Customer Service Department at (203) 377-8282. Have your purchase order number available before calling Newport. The Customer Service Representative will give you a Return Material Authorization number (RMA). Having an RMA will shorten the time required for repair, because it ensures that your equipment will be properly processed. Write the RMA on the returned equipment's box. Equipment returned without a RMA may be rejected by the Newport Receiving Department. Equipment returned under warranty will be returned with no charge for the repair or shipping. Newport will notify you of any repairs not covered by the warranty, with the cost of the repair, before starting the work.

Please return equipment in the original (or equivalent) packaging. You will be responsible for damage incurred from inadequate packaging, if the original packaging is not used.

Include the cables, connector caps and antistatic materials sent and/or used with the equipment, so that Newport can verify correct operation of these accessories.